# MICRODIRECT™ 1080p HD

HANDHELD DIGITAL MICROSCOPE

ITALIAN

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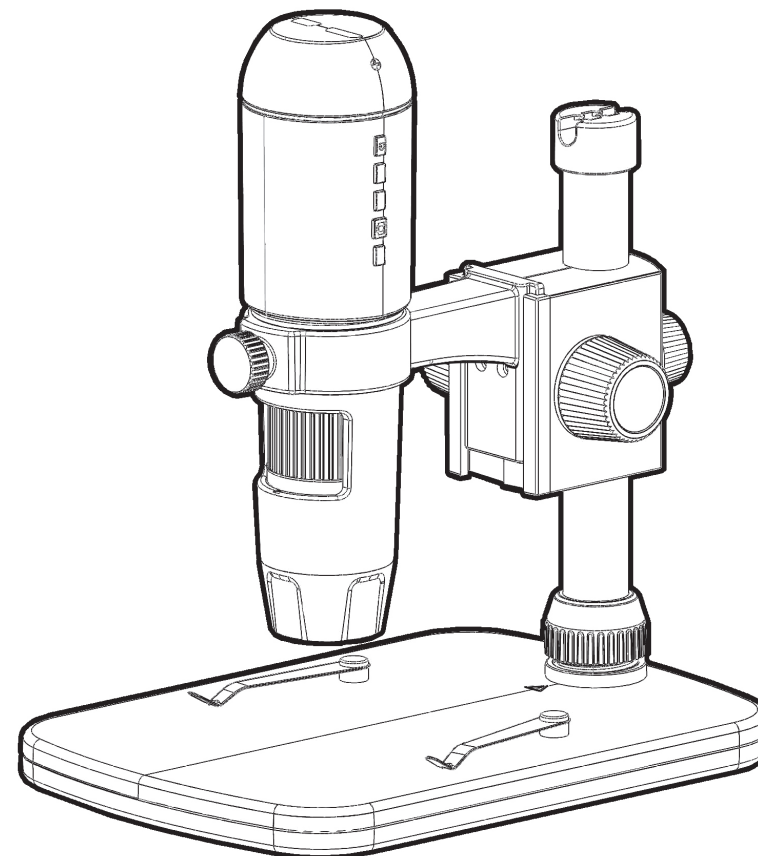
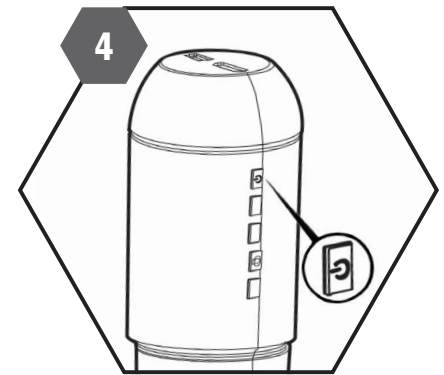
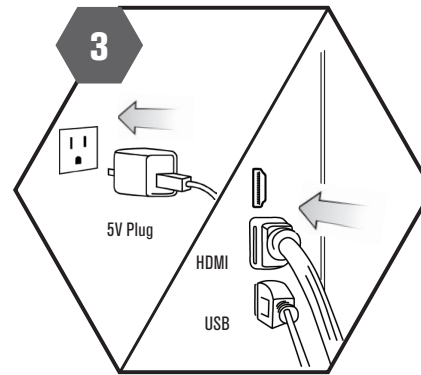
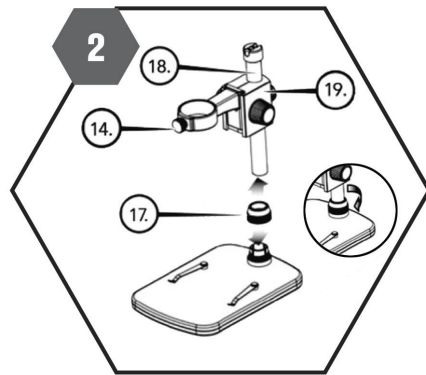
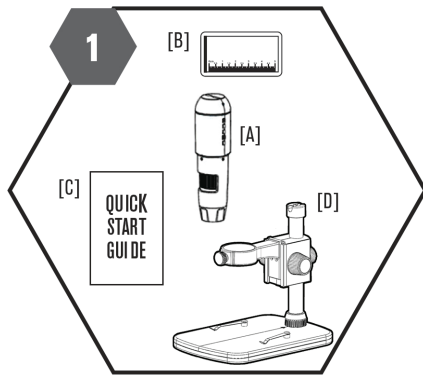
ENGLISH

### 1. WHAT IS IN THE BOX

- [A] 1 x MicroDirect\* 1080p HD Microscope
- [B] 1 x Calibration Ruler
- [C] 1 x Quick Start Guide
- [D] 1 x Portable Stand
- [E] 1 x HDMI Cable
- [F] 1 x USB Cable

### 2. SETUP THE MICROSCOPE AND STAND

- Carefully remove all components from the box.
- The microscope holder assembly (19) should be connected to the support shaft (18) in the box. If it is not, connect it now.
- Attach complete assembly (17, 18, 19) to the base threads and tighten.
- Insert the microscope into the microscope holder and tighten the microscope holder lock nut (14) in place.

**NOTE:** The microscope can be used in handheld mode as well without the stand.

### 3. CONNECT HDMI AND USB TO MONITOR OR PROJECTOR

- Connect the included HDMI cable to the microscope and to the open HDMI port on the monitor or projector.
- Plug the USB cable into the microscope and then into an open USB 2.0 port on your monitor or projector (if available).

**NOTE:** If the monitor or projector you are connecting to does not have a standard 2.0 USB plug- with power, you must use an alternative power source. Suggestions are:

- Plug into any 5v-1A to 2.1A plug and plug into 120V outlet.
- Use a 5V power bank, like Celestron Elements series, with an open USB port.

[www.celestron.com/elements](http://www.celestron.com/elements)

### 4. TURN MICRODIRECT® 1080P HD MICROSCOPE ON

- Press the ON/OFF button on the microscope.

For the complete instruction manual go to this product's page at [www.celestron.com](http://www.celestron.com) and click on the SUPPORT section.

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